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Cat number: KGY014

Store at : -20°C for 12 months

For Research Use Only (科研专用)

Catalog Number:	KGY014
Content:	50mg Plasmocin™ Treatment, provided in 2 x 1 ml tubes at a concentration of 25 mg/ml.
Storage Condition:	Store at -20°C
Shelf Life:	Stable for up to 12 months at -20°C. Before storing at -20°C, prepare aliquots to avoid repeated freeze-thaw cycles.
Note:	Presence of crystals does not alter properties of the product. Vortex the tube until the crystals disappear.
Method:	<p>Treatment of Mycoplasma Infect Cell Culture:</p> <p>Typically, Plasmocin Treatment is used for 2 weeks at 25 µg/ml which represents a 1:1000 dilution of the 25mg/ml stock solution (100µl Plasmocin Treatment in 100ml culture medium).Plasmocin™ Treatment. can be used in combination with penicillin and streptomycin.</p> <ol style="list-style-type: none">1.Split an actively dividing culture of cells into medium containing 25 µg/ml of Plasmocin™ Treatment.2. Remove and replace with fresh Plasmocin™ Treatment containing medium every 3-4 days for 2 weeks. <p><i>Note: If 25µg/ml of Plasmocin is toxic for your cells, reduce the concentration of Plasmocin™ to 12.5 µg/ml and treat for an additional week.</i></p> <ol style="list-style-type: none">3. Test for the presence of mycoplasma by a conventional PCR method or using a cell based assay. If mycoplasmas are still present following treatment, see troubleshooting overleaf. <p><i>Note: InvivoGen provides Plasmotest a simple, rapid and reliable assay for the visual detection of mycoplasma contamination. This is the first assay to utilize cells to signal the presence of mycoplasma.</i></p> <ol style="list-style-type: none">4. For the maintenance of a mycoplasma free culture, use Plasmocin™ Prophylactic at a concentration of 5 µg/ml.

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